



## Acute 3D hiPSC-CM Contractility Tool

### PROTOCOL:

#### 1. Preparation of plates and media

- 1.1. In a sterile tissue culture hood, prepare a sterile 6-well plate by transferring 2 mL of 0.1% gelatin (**Table of Materials**) to each well. Place the lid on the 6-well plate and allow the coated plate to incubate at 37 °C for a minimum of 1 h.
- 1.2. One day before seeding hiPSC-CMs and primary human ventricular cardiac fibroblasts on ECT (Engineered cardiac tissue) molds to form 3D tissues, thaw hydrogel (**Table of Materials**) aliquot in the refrigerator, on ice.
- 1.3. Prepare cardiomyocyte medium (**Table of Materials**) according to the manufacturer's instructions.
- 1.4. Prepare cardiac fibroblast medium (**Table of Materials**) according to the manufacturer's instructions.

#### 2. Seeding of cryopreserved hiPSC-CMs

- 2.1. Two days before seeding hiPSC-CMs on ECT molds, pre-plate hiPSC-CMs on 0.1% gelatin-coated, sterile 6-well plates. Thaw hiPSC-CMs using standard thawing protocol <sup>1,2</sup>.
- 2.2. Then, plate 1,500,000 total hiPSC-CMs (**Table of Materials**) per well according to the manufacturer's instructions <sup>3</sup>.
- 2.3. Culture hiPSC-CMs in standard cardiomyocyte media for 2 days to allow hiPSC-CMs to recover from cryopreservation at 37 °C, 5% CO<sub>2</sub>. Refresh the supernatant with 100% cardiomyocyte medium every 48 h.

#### 3. Seeding of cryopreserved cardiac fibroblasts

- 3.1. Two days before seeding cardiac fibroblasts on ECT molds, pre-plate cardiac fibroblasts on 0.1% gelatin-coated, sterile 6-well plates. Thaw cardiac fibroblasts using standard thawing protocol <sup>1,2</sup>.
- 3.2. Then, plate 250,000 total cardiac fibroblasts (**Table of Materials**) per well according to the manufacturer's instructions <sup>3</sup>.
- 3.3. Culture cardiac fibroblasts in standard cardiac fibroblast medium for 2 days at 37 °C, 5% CO<sub>2</sub>. Refresh the supernatant with 100% cardiac fibroblast medium every 48 h.

#### 4. Dissociation and counting pre-plated hiPSC-CMs and cardiac fibroblast.

- 4.1. Check the status of hiPSC-CMs before dissociation. Evaluate hiPSC-CM health ensuring viability and stable beating.

NOTE: The purity of the hiPSC-CM population is important (e.g., >90 % Cardiac Troponin T) <sup>4</sup>. A cardiomyocyte selection method (e.g., metabolic selection or sorting) is recommended to reduce



disruption by non-cardiomyocyte cells <sup>2,3</sup>.

- 4.2. Wash hiPSC-CMs 2x with 4 mL per well of D-PBS without CaCl<sub>2</sub> or MgCl<sub>2</sub> (**Table of Materials**). Aspirate D-PBS and add 1 mL of room temperature dissociation reagent to each well then incubate for 15 min at 37 °C<sup>5</sup>.
- 4.3. Add 10 mL of cardiomyocyte medium to a sterile 15 mL conical tube.
- 4.4. Dissociate hiPSC-CMs from the 6-well plate with a 1,000 µL pipette (**Figure 1B**). Add the cell suspension to the 15 mL conical tube <sup>3</sup>.
- 4.5. Rinse the well with 1 mL of fresh cardiomyocyte medium to collect any residual hiPSC-CMs and add to the 15 mL conical tube. Bring the final volume of the conical tube to 15 mL.
- 4.6. Centrifuge for 5 min (200 × *g*). Remove the supernatant up to the 1 mL mark. Resuspend the cells in cardiomyocyte medium to a final volume of 5 mL.
  - 4.6.1. Count hiPSC-CMs with a manual or automated cell counter.
- 4.7. Incubate the hiPSC-CM suspension at room temperature while the cardiac fibroblast are dissociated (30 min maximum).
- 4.8. Dissociate and count cardiac fibroblast as above.

## 5. Generation of Engineered Cardiac Tissues<sup>6</sup>

- 5.1. Combine hiPSC-CMs and cardiac fibroblasts at a 10:1 ratio (i.e., 100,000 hiPSC-CM:10,000 cardiac fibroblast)<sup>5,6</sup>.
- 5.2. Combine cell mixture and Extracellular Matrix hydrogel components (**Table of Materials**).
- 5.3. Seed cell-hydrogel suspension into each well of ECT molds.
- 5.4. Culture tissues for 7 days to enable remodeling and compaction at 37 °C, 5% CO<sub>2</sub> refresh medium every 48 h.
- 5.5. Expose tissues to 7-week electrical condition stimulation protocol<sup>5,6</sup>.

NOTE: Electrical stimulation protocol of weekly 1 Hz increase in frequency is recommended for ventricular functional maturation until positive force-frequency response is observed<sup>5,6</sup>.

## 6. Contraction recording and analysis<sup>7</sup>

- 6.1. Prepare CCM assay medium, Tyrode's solution containing (in mmol/L): CaCl<sub>2</sub> 1.8, NaCl 134, KCl 5.4, MgCl<sub>2</sub> 1, glucose 10, and HEPES 10, pH adjusted to 7.4 with NaOH, and equilibrate to 37 °C in a water bath.
- 6.2. Place tissue in tissue bath (1 cm × 5 cm, 600 µl) recording chamber.



- 6.3. Perfuse 37 °C CCM assay medium at 600  $\mu$ L per chamber at a flow rate of 4 ml/min.
- 6.4. Cut and attached one end of the tissue to a force transducer (**Table of Materials**)<sup>7</sup>.
- 6.5. Immobilize the other end of the tissue using stainless steel wires attached to a micromanipulator<sup>7</sup>.
- 6.6. Connect a silicon-based strain gauge with two piezoresistive elements (**Table of Materials**) to the amplifier to convert resistance change to voltage signals. Convert voltage to force measurements.
- NOTE: Prior to experimentation the strain gauge was pre-calibrated with known weights and a conversion factor of 104.4  $\mu$ N/mV was established<sup>7</sup>.
- 6.7. Use camera (**Table of Materials**) to capture contraction videos and force transducer to measure contractility. Field stimulate the tissue with a commercial pulse generator (**Table of Materials**) to electrically pace the tissue. Pace tissue at 1.5x threshold at 1 Hz with baseline pulse parameters (e.g., monophasic square wave pacing pulses with a 2 ms stimulus pulse duration ( $\sim$ 10 V/cm)<sup>7,8</sup>.
- 6.8. Record the baseline, pacing only (i.e., before CCM) contraction for a minimum of 2 min<sup>8,9</sup>.
- 6.9. Then, stimulate the tissue with an experimental electrical signal. To follow this protocol, use the standard CCM stimulation parameters: two symmetrical biphasic pulses of **5.14 ms phase duration (20.56 ms total duration)**,  **$\sim$ 28 V/cm (phase amplitude)**, **zero interphase interval** and **30 ms delay** (i.e., time from the end of the pacing pulse and the beginning of the CCM pulse)<sup>10,11</sup> and record CCM-induced contraction for a minimum of 2 min<sup>7</sup>.
- 6.10. Turn off the CCM signal and stimulate with baseline pacing pulse and record contraction of the recovery period (i.e., after CCM) for a minimum of 2 min.
- 6.11. Enable contractile properties to return to baseline then use contractility tool to evaluate various experimental electrical signals.
- 6.12. Use standard contraction software to analyze contraction videos and conversation factor to quantify contraction force (e.g., **contraction/force amplitude, contraction slope, relaxation slope, time to peak, time to baseline 90%**, and **contraction duration 50%**)<sup>4,7,8,12,13</sup>.

#### Table of Materials

Name of Material/ Equipment	Company	Catalog Number	Comments/Description
0.1% Gelatin	STEMCELL Technologies	7903	Pre-plating Culture Substrate
6-well Plate	Thermofisher	140675	hiPSC-CM Culture, Plastic
Calcium Chloride dihydrate (CaCl <sub>2</sub> )	Fisher Scientific	c70-500	Tyrode's solution
Conical tube 15 mL	Corning	352099	Dissociation
Digital CMOS Camera	Hamamatsu	C11440-42U30	Contraction Video Recording



D-PBS	Life Technologies	14190-144	Cell Wash
Glucose	Sigma-Aldrich	G8270-1kg	Tyrode's solution
Hemocytometer	Fisher Scientific	22-600-107	Cell Counting
HEPES	Sigma-Aldrich	H3375	Tyrode's solution
iCell Cardiomyocytes Plating Medium	Fujifilm Cellular Dynamic, Inc.	M1001	hiPSC-CM Plating Media
iCell Cardiomyocytes <sup>2</sup> , 01434	Fujifilm Cellular Dynamic, Inc.	R1017	hiPSC-CMs
Incubator (37 °C, 5% CO <sub>2</sub> )	Thermofisher	50116047	Maintain cells and tissues
Inverted Microscope	Olympus	IX73	Imaging ECTs
Magnesium Chloride hexahydrate (MgCl <sub>2</sub> )	Fisher Scientific	m33-500	Tyrode's solution
Matrigel Basement Membrane Matrix	Corning	354230	ECM Component
Microcentrifuge tubes 1.5 ml	Fisher Scientific	05-408-129	Substrate Aliquot
Model 3800 Multichannel Power Stimulator	AM-Systems	Model 3800	Pulse Generator
Pen-Strep	Invitrogen	15140-122	Cardiomyocyte Media
Pipette L-20	Rainin	17014392	Pipette
Pipette P1000	Fisher Scientific	F123602G	Pipette
Pipette tips, 1000 ul	Fisher Scientific	02-707-509	Pipette
Pipette tips, 20 ul	Rainin	GPS-L10S	Pipette
Potassium Chloride (KCl)	Fisher Scientific	P330-500	Tyrode's solution
Sodium Chloride (NaCl)	Fisher Scientific	s641-212	Tyrode's solution
Sodium Hydroxide (NaOH)	Sigma-Aldrich	221465	Tyrode's solution
Stimulation Electrodes	--	--	Pacing and CCM Stimulation
Trypan Blue Stain	Life Technologies	T10282	Cell Counting
TrypLE Express	Life Technologies	12605-010	Cell Dissociation
Fibrin	Sigma-Aldrich	9001-31-4	ECM Component
Collagenase type 2	Worthington	4176	ECM Component
Multi-Purpose Sensor Element	Kronex Technologies	AE801	Silicon-based strain gauge/force transducer
NHCF-V – Human Ventricular Cardiac Fibroblasts	Lonza Bioscience	CC-2904	Cardiac fibroblasts
FGM™-3 Cardiac Fibroblast Growth Medium-3 BulletKit	Lonza Bioscience	CC-4526	Cardiac fibroblast medium



## REFERENCES:

- 1 Feaster, T. K. *Implementation of human-induced pluripotent stem cell-derived cardiomyocyte to model excitation-contraction coupling in health and disease* Doctor of Philosophy thesis, Vanderbilt University Medical Center, (2015).
- 2 Ma, J. *et al.* High purity human-induced pluripotent stem cell-derived cardiomyocytes: electrophysiological properties of action potentials and ionic currents. *Am J Physiol Heart Circ Physiol.* **301** (5), H2006-2017, (2011).
- 3 Burridge, P. W., Holmstrom, A. & Wu, J. C. Chemically Defined Culture and Cardiomyocyte Differentiation of Human Pluripotent Stem Cells. *Curr Protoc Hum Genet.* **87** 21 23 21-21 23 15, (2015).
- 4 Feaster, T. K. *et al.* Matrigel Mattress: A Method for the Generation of Single Contracting Human-Induced Pluripotent Stem Cell-Derived Cardiomyocytes. *Circ Res.* 10.1161/CIRCRESAHA.115.307580, (2015).
- 5 Zhao, Y. *et al.* A Platform for Generation of Chamber-Specific Cardiac Tissues and Disease Modeling. *Cell.* **176** (4), 913-927.e918, (2019).
- 6 Sun, X. & Nunes, S. S. Maturation of Human Stem Cell-derived Cardiomyocytes in Biowires Using Electrical Stimulation. *J Vis Exp.* 10.3791/55373 (123), (2017).
- 7 Feaster TK, F. N., Pallotta I, Narkar A, Casciola M, Graziano MP, Aschar-Sobbi R and Blinova K. Acute effects of cardiac contractility modulation stimulation in conventional 2D and 3D human induced pluripotent stem cell-derived cardiomyocyte models. *Front. Physiol.* **13:1023563**, (2022).
- 8 Feaster, T. K., Casciola, M., Narkar, A. & Blinova, K. Acute effects of cardiac contractility modulation on human induced pluripotent stem cell-derived cardiomyocytes. *Physiol Rep.* **9** (21), e15085, (2021).
- 9 Narkar, A., Willard, J. M. & Blinova, K. Chronic Cardiotoxicity Assays Using Human Induced Pluripotent Stem Cell-Derived Cardiomyocytes (hiPSC-CMs). *Int J Mol Sci.* **23** (6), (2022).
- 10 ImpulseDynamics. *OPTIMIZER® Smart Implantable Pulse Generator INSTRUCTIONS FOR USE*, <<https://impulse-dynamics.com/wp-content/uploads/2020/05/13-290-008-01-US-Rev-01-OPT-Smart-IPG-IFU.pdf>> (2018).
- 11 ImpulseDynamics. *OPTIMIZER™ Smart Mini Implantable Pulse Generator INSTRUCTIONS FOR USE*, <<https://impulse-dynamics.com/global/wp-content/uploads/sites/2/2021/02/13-290-011-EU-Rev-00-OPTIMIZER-Smart-Mini-IPG-IFU-EU.pdf>> (2019).
- 12 Sala, L. *et al.* MUSCLEMOTION: A Versatile Open Software Tool to Quantify Cardiomyocyte and Cardiac Muscle Contraction In Vitro and In Vivo. *Circ Res.* **122** (3), e5-e16, (2018).
- 13 Grune, T., Ott, C., Haseli, S., Hohn, A. & Jung, T. The "MYOCYTER" - Convert cellular and cardiac contractions into numbers with ImageJ. *Sci Rep.* **9** (1), 15112, (2019).